Comparison of Two Analytical Chemistry Techniques

A Lab Experience for High School
Chemistry Students

Abigail Weaver Clay High School aweaver@sbcsc.k12.in.us



Teaching Goals

- Students would be introduced to acid base back titrations as a tool to measure concentration of diethlycarbamazine (DEC) citrate.
- Students would be introduced to a method of determining accuracy of a procedure.

• Students would learn a technique to compare two sets of data.



The Role of DEC in Fighting a Tropical Disease.

Man with limb swollen due to effects of LF. Carter Center Photo: E. Staub

- Lymphatic Filariasis (LF) affects 120 million people with more than a billion at risk. Over 40 million people are permanently disabled by this disease making it a leading cause of disability in the world. Consuming salt fortified with diethylcarbamazine (DEC) has been found to be a safe and effective method of prevention for LF¹.
- This titration was developed at the University of Notre Dame and is used for quality control of DEC citrate fortified salt in Haiti
- 1. Fifth Meeting of the Global Alliance to Eliminate Lymphatic Filariasis. Rep. The Global Alliance to Eliminate Lymphatic Filariasis, Apr. 2008. Web. 24 July 2009. http://www.filariasis.org/documents/GAELF5reportEnglish.pdf

Student Activities



Bon Sel salt contains DEC citrate to prevent lymphatic filariasis

- Produce a standard curve by titration of known samples.
- Measure DEC citrate concentration using the acid base back titration method
- Comparison of measured concentrations vs. actual concentrations
- Comparison of DEC citrate concentrations measured by titration method vs. HPLC method





A student titrating DEC citrate.

The Titration Procedure

- An acid-base back titration is one method used to quantify DEC citrate
- The titration method uses the following sample:


 10.00mL solution of DEC citrate in 10% saline
 10.00mL of 0.01M sodium hydroxide
 2 drops phenolphthalein
- The sample is titrated with 0.004M hydrochloric acid to a clear endpoint.



Student Lab Work

- 1) Using samples of known DEC citrate concentration students can produce a standard curve.
- 2) Using the standard curve students can determine the concentration of DEC citrate in titrated unknowns.

DEC Citrate Standard Curve

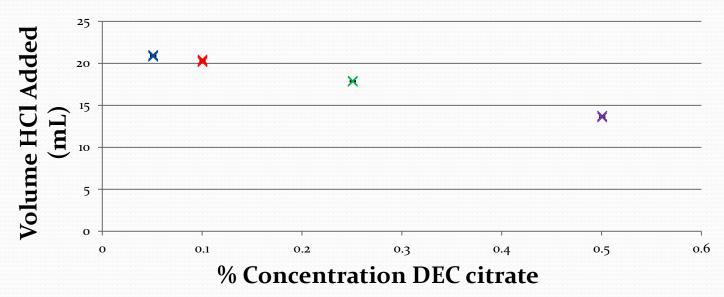




Accuracy of DEC Citrate Titration Data analysis

*Students can graph their titration concentrations versus the actual concentration to estimate accuracy.


*A slope of 1.00 indicates accurate measurements.

Actual vs. Experimental % DEC Citrate Concentration

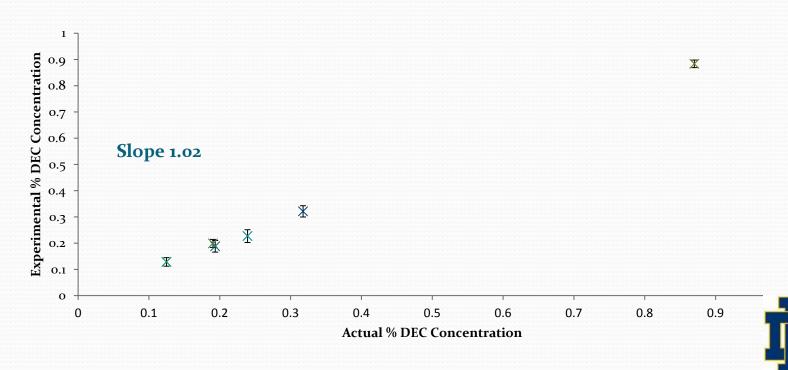

Comparison of Two Methods

• Students can compare their results to measurements acquired by another technique.

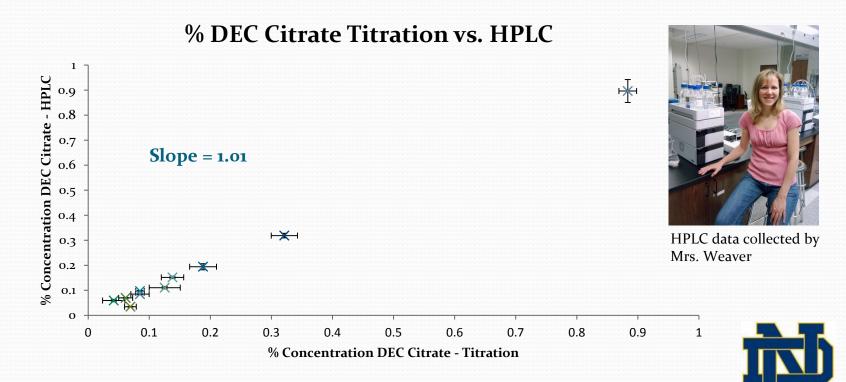

Conclusions



A Haitian chemist testing DEC fortified salt Photo: Sarah Miller, University of Notre Dame

- Students gain experience with an acid base back titration
- Students learn methods to evaluate an analytical procedure
- Students gain experience graphing with EXCEL
- Students learn real world application of a titration
- Concentrations of compounds used pose no health or safety concerns to students

Thanks to The University of Notre Dame Research Experience for Teachers Program, The Notre Dame Haiti Program, The Lieberman Group, Doug Miller, Shannon Huey, Béatrice Blanc and Steve Wietstock

